PCE Italia s.r.l.
Via Pesciatina
878 / B-Interno 6
Loc. Gragnano - Capannori
55010 Lucca (Italia)
Tel.: +39 0583 975 114
Fax: +39 0583 974 824
info@pce-italia.it
www.pce-italia.it

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OPERATION
Place scale on a horizontal flat surface.
1. Place scale on a horizontal flat surface, press 'ON/OFF' key.
2. Wait until '0.0' is displayed.
3. Put the objects on the weighing platform.
4. Press 'CAL MODE' key, you can switch among the weighing units.

TARE WEIGHING
1. Turn off scale as described above.
2. Place the same item on the platform.
3. Press 'TARE' key and wait until '0.0' is displayed.
4. Add the 'net-weight/item'.

CALIBRATION
The gravity change from place to place in the world, you may need to calibrate your scale to adjust it to the local gravity by pressing 'ON/OFF' key.
1. Turn on the scale by pressing 'ON/OFF' key.
2. Press 'CAL MODE' key and hold it until LCD display 'CAL' and followed by a number which is the weight you need to put on the platform.
3. Check 'MODE' key again, the 'CAL' will flash and followed by a number which is the weight of the required standard list, weight on the platform and wait 3-4 seconds, LCD will show 'pass'.

Special Calibration
The gravity change from place to place in the world, you may need to calibrate your scale to adjust it to the local gravity by pressing 'ON/OFF' key.
1. Turn on the scale by pressing 'ON/OFF' key.
2. Press 'CAL MODE' key and hold it until LCD display 'CAL' and followed by a number which is the weight you need to put on the platform.
3. Check 'MODE' key again, the 'CAL' will flash and followed by a number which is the weight of the required standard list, weight on the platform and wait 3-4 seconds, LCD will show 'pass'.

Special Calibration
1. Turn off the scale.
2. Press both 'CAL' and 'TARE' key until 'CAL' is displayed on screen.
3. Release 'CAL' and 'TARE' key, the internal number will be displayed on LCD.
4. Press 'CAL' key again, the display will show '2000' and then a number.
5. Put the required standard list weight on the inside platform, wait 3-4 seconds.
6. Press 'CAL' key and press 'ON/OFF' key to turn off the scale.
7. Remove the weight and press 'ON/OFF' key to calibrate the scale by the above steps.
If the display shows 'FAIL', please re-calibrate the scale by the above steps.

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